

## Adam Equipment

## PT PLATFORM SET-UP

Thank you for purchasing your Adam Equipment AE PT Platforms.

The Scale is a single base with one loadcell cable connection.

- 1) To remove the scale from the packaging, attatch the eye bolts supplied and lift the platform using a suitable lifting device, i.e. fork truck or gantry.
- 2) Fit the 4 feet supplied to the underneath of the platform.
- 3) Place the platform on the ground ensuring that the platform is level using the 4 feet to adjust and using the built-in spirit level make sure that the platform is stable and does not rock.
- 4) Take care to make sure that the cable is not rapped or pinched and is not causing a trip hazard.
- 5) Connect the indicator end of the cable to your indicator and follow the indicator instructions.