




SCIOLOGEX

SCI-Stir

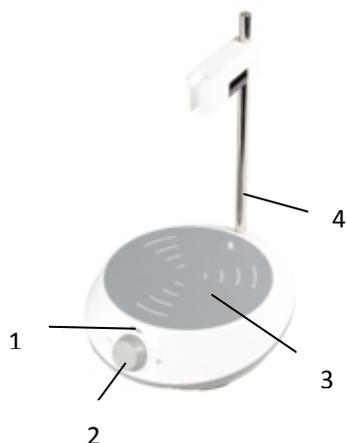

1. POWER LED
2. SPEED ADJUSTMENT KNOB
3. WORK PLATFORM
4. SENSOR HOLDER

SCIOLOGEX, LLC, 1275

1275 Cromwell Ave.

Suite C6, Rocky Hill, CT 06067 USA

Tel: +1(860) 436-9221

Fax: +1(860) 436-9745

info@scilogex.com

www.scilogex.com

Item	Parameter
Voltage [VAC]	100-240
Frequency [Hz]	50/60
Power Input [W]	1.2W
Adjustment	Scale
Speed Range [rpm]	300—2000
Max. Capacity [L]	1.5
Max. Stir Bar	35mm
Work Platform [mm]	Ø120mm
Work Platform Material	PET
Casing Material	ABS
Dimension D×W×H [mm]	145×150×50
Weight [kg]	0.3
Permissible Temperature [°C]	5-40
Permissible Humidity	80%
Protection Class	IP21

Safety Precautions

Please read carefully before operating this instrument.

- Ensure that only trained staffs use the instrument.
- Never use the instrument in any manner not specified in these instructions.
- If any abnormal condition occurs during operation, please power-off immediately and contact our service center.
- **Caution** – Effects of the magnetic field have to be taken into consideration (e.g. data storage media, cardiac pacemakers).
- Please avoid splashes/spills.

Operation

Place the EcoStir on a level surface; confirm the EcoStir is working under the rated voltage. Rotate the speed adjustment knob to control the EcoStir stirring speed, clockwise to accelerate, anti-clockwise to decelerate.

EcoStir's speed can be adjusted from 300rpm to 2000rpm.

Warranty

This product is guaranteed for two years.